

HOW TO SWITCH FROM DAYBOOK 2 (2018-2019) TO DAYBOOK 3 (2020)

Daybook 3 is a new software released in April 2020.

Many improvements have been made in the features and algorithms precision.

The results of Daybook 2 cannot be transferred directly into Daybook 3.

You will have to recreate in Daybook 3 (within the Data Manager module) your system, your acquisition profile and repopulate your timeline.

1- How to install Daybook 3

You do not need to uninstall Daybook 2.

1) Download Daybook 3 from our website www.argolight.com

2) **Make sure to fully close any session of Daybook 2.**

("Mongodb" and "node" process should not be running within the task manager)



Daybook 2 and Daybook 3 should not be launched on the same computer at the same time

3) Launch the install.

4) Open Daybook 3

5) On the top left corner, use the "Register" button to enter your license code.

For any question regarding license purchase or renewal, please write to

contact@argolight.com

2- How to input your data in Daybook 3

A. Easily create your microscope:

Make sure to close any session of Daybook 2 before launching Daybook 3.

Use the "Import from image" button within the Data Manager module to pre-fill the fields directly from a proprietary image file. The documentation within Daybook 3 (Data Manager module) gives a detailed explanation on this feature.

B. Easily create your acquisition profile:

Use the "Autofill the form with image data" button to pre-fill the fields directly: simply verify the filled-in fields and/or complete the blank ones.

C. Easily re-analyze your images:

Use the "Batch Analysis" button within the Analysis module to run several analyses in one go on a batch of raw images. The documentation within Daybook 3 (Analysis module) gives a detailed explanation on this feature.

A trouble to install Daybook 3? customer@argolight.com

More info about Daybook 3 price and options? contact@argolight.com